

DeepPIV on MiniROV Deployment #10 Checklist

v 1.2.0

11Nov15

K. Katija

Dive Objective: Deep PIV flow & particle measurements around
Bathochordus stigmus & B. minutus houses. Also looking for other
relatively organisms for 3D reconstruction proof-of-concept.
Fluorescent dye will be used to elucidate bulk flow properties.
(MiniROV dive #62)

Deployment Date: 12-01-15

Recovery Date: _____

Person performing topside checklist: KAKANI KATIJA

Pre-deployment:

1. Ensure that comm and power systems on MiniROV are functioning normally.
2. Align the laser and camera
 - N/A [a. Attach protective shroud and cloth to MiniROV's body
 - i. Zero digital level based on orientation of MiniROV sitting on deck
 - ii. Verify alignment of shroud using a digital level
 - iii. Record digital level reading on shroud: _____
 - ✓ b. Enable Tilt and Laser controls in LabVIEW
 - ✓ c. Zero the Science Camera
 - i. Click Zero button in Tilt dialog tab
 - ii. Click Home button in Tilt dialog tab
 - iii. Record digital level reading on camera: _____
 - ✓ d. Remove laser housing protective cover, and put on protective goggles
 - ✓ e. Power up the laser
 - i. Rotate the dial to 1.5 V in Laser dialog tab
 - ii. Record output of ADC: _____
 - N/A [f. Verify that laser sheet is perpendicular to Science camera
 - i. Measure angle of laser sheet on shroud using a digital level
 - ii. Record digital level reading of laser sheet: _____
 - ✓ g. Power down the laser
 - h. Replace laser housing protective cover
- N/A [3. Remove protective shroud from MiniROV.
4. Secure protective shroud on deck of ship.
5. Run pre-dive test program and verify correct operation of the following systems:
 - ✓ a. MiniROV LED lights
 - ✓ b. Science camera (adjust zoom, focus, shutter speed, and IR filter settings)
 - ✓ c. AJA HD recorder (record, stop, playback)
 - ✓ d. Dye injector/vortex generator (disconnect tubing if dye injector being used, move piston, reset piston, set different increments)
 - ✓ e. Sizing lasers (measure separation at multiple distances)
- ✓ 6. Sync all components (including Science camera) with ship's NTP

7. Start MiniROV dive log
8. Check that all cable connectors and dummies are installed and tight.
9. Check that all dye/vortex generator tubing is installed and tight.
10. Open all dye valves and connectors.
11. Remove laser housing protective cover.
12. Check that the MiniROV tool sled is secure.
13. Check for tools left on and in MiniROV frame.
14. Remove MiniROV hold-down straps.

Deployment:

1. Start AJA recorder
2. Deploy MiniROV, noting exact time and position of launch.
 - a. MiniROV launch time: 15:38:00 PDT
 - b. Latitude: 36° 44.76 deg
 - c. Longitude: 122° 6.297 deg
3. Record time it takes for MiniROV to submerge: 15:39:36
4. Dive plan: refer to _____

Recovery:

5. Check that the laser is powered down.
6. Change the Science camera settings (focus and others) to auto and zoom to wide.
7. Complete MiniROV dive log
8. Stop recording video with Aja
9. Recover MiniROV
10. Secure MiniROV and DeepPIV on the ship deck.
11. Record the time that MiniROV is on deck: 18:35 PDT
12. Close all dye injector/vortex generator valves.
13. Replace laser housing protective cover.