

DeepPIV on 04.05.17 PAULA CARSON Deployment # 57 Checklist
v 5.0.0 22 August 16
K. Katija

Dive Objective: MICROPLASTICS

- plastic injection w/ salt chain
- DeepPIV attempt w/ batho

Deployment Date: 04.05.17
Recovery Date: 04.05.17
Person performing topside checklist: KARSON

Pre-deployment alignment:

1. Verify that DeepPIV laptop and peripherals are connected (Ethernet, usb, mouse, monitor, AJA/Shogun recorder)
2. Power up DeepPIV
3. Open Silex USB Controller and connect
4. Open Sony Remote Camera Controller and modify camera settings
5. Verify laser and camera alignment
 - a. Place protective shroud around DeepPIV instrumentation
 - b. Remove protective covers on laser and camera housings
 - c. Check that housing markings on laser and camera housing are aligned
 - d. Power up the laser
 - i. Rotate the dial to 2 V in Laser dialog tab
 - ii. Record output of ADC: _____
 - iii. Measure laser power: _____
 - e. Using digital level, compare laser sheet alignment with camera housing glass port
 - i. Record digital level reading on laser sheet: _____
 - ii. Record digital level reading on camera housing: _____
 - f. Power down the laser
 - g. Replace protective covers on laser and camera housings
 - h. Remove protective shroud from DeepPIV instrumentation
 - i. Close Remote Camera Controller software, disconnect Silex USB connection
6. Power down DeepPIV

Pre-deployment:

- ✓ 1. Verify that DeepPIV laptop and peripherals are connected (Ethernet, USB, mouse, monitor, AJA/Shogun recorder)
- ✓ 2. Power up DeepPIV
- ✓ 3. Run pre-dive test program and verify correct operation of the following systems:
 - ✓ a. DeepPIV camera (ISO, shutter speed, f-stop, record)
 - i. Open Silex USB Controller and connect
 - ✓ ii. Open Sony Remote Camera Controller
 - b. DeepPIV laser

- i. Power up the laser, rotate the dial to 2 V
 - ii. Record output ADC: 2014
 - iii. Measure laser power: N/A
- ✓ c. AJA/Shogun HD recorder (record, stop, playback)
- ✓ d. Dye injector/vortex generator (disconnect tubing if dye injector being used, move piston, reset piston, set different increments)
 - i. Note COM port on ROV: _____
 - ii. Note COM port on Serial-to-USB converter (either 1/2): _____
 - iii. Open ELMO Controller software
 - iv. Direct connect to motor
 - v. Set baud rate to 19200
 - vi. Set COM port (either 3/4): _____
 - vii. Connect
 - viii. Use Velocity Mode, hit stop to prevent dye leakage
 - ix. Set speed to 100 rpm and hit go, stop; reverse direction, go, stop
- e. Translational stage
 - i. ~~Ensure that Ethernet cable is connected~~
 - ii. ~~Open ELMO Controller software~~
 - iii. ~~Direct connect to motor~~
 - iv. ~~Set baud rate to 19200~~
 - v. ~~Set COM port (either 1/2): _____~~
 - vi. ~~Connect~~
 - vii. ~~Use Velocity Mode, hit stop to prevent screw from turning~~
 - viii. ~~Set speed to 500 RPM and hit go, stop; reverse direction, go, stop~~
- f. Sizing lasers
- ✓ 4. Sync all components (including Science and DeepPIV camera) with ship's NTP and record video on both cameras
- ✓ 5. Check that all cable connectors and dummies are installed and tight
- ✓ 6. Check that all dye/vortex generator tubing is installed and tight
- ✓ 7. Open all dye valves and connectors
- ✓ 8. Remove protective covers on laser and camera housings

Deployment:

1. Deploy ROV, noting exact time and position of launch.
 - a. ROV launch time: _____ PDT
 - b. Latitude: 30°11.9022 deg
 - c. Longitude: 122°3.0900 deg
2. Record time it takes for ROV to submerge: _____
3. Dive plan: refer to _____
4. Start AJA/Shogun recorder when event of interest occurs; stop when finished

Recovery:

1. Record the time that ROV is on deck: _____ PDT
2. Power down DeepPIV
3. Close all dye injector/vortex generator valves
4. Rinse instrument housings and attachments with fresh water
5. Replace protective covers for laser and camera housings