

User's Manual ATH 31

Introduction

■ Introduction to the ATH 31 pump and ACT 201 H controller	■ A 10
■ The hybrid turbomolecular pump operating principle	■ A 20
■ The different versions	■ A 30
■ The ACT 201 H controller	■ A 40
■ Integrable ACT 201 H "mini Board" controller . .	■ A 50
■ The accessories	■ A 60
■ The technical characteristics	■ A 70
■ Hybrid pump dimensions	■ A 80
■ Integrable ACT 201 H "mini Board" controller technical characteristics	■ A 91

Start-up

■ Safety instructions	■ B 10
■ Pump connection to an installation	■ B 20
■ Water cooling circuit connection	■ B 30
■ Nitrogen purge connection for corrosive version .	■ B 40
■ Venting valve connection	■ B 50
■ Air cooling kit connection	■ B 60
■ ACT 201 H electrical connections	■ B 70
■ «Remote Control» connector wiring (ACT 201 H box version)	■ B 80
■ ACT 201 H "mini Board" power supply and remote control wiring	■ B 81
■ RS 232 or RS 485 serial link wiring (ACT 201 H box version)	■ B 90
■ RS 232 or RS 485 serial link wiring (ACT 201 H "mini Board")	■ B 91
■ Detailed description of RS commands	■ B 100

User's Manual ATH 31

Start-up (continued)	■ The controller start-up	■ B 110
	■ Configuring the controller for the application . . .	■ B 120
	■ First pump start-up	■ B 130
Operation	■ Hybrid-turbo pump in a pumping application . .	■ C 10
	■ Controlling the pump using the controller front panel.	■ C 20
	■ «Remote Control» connector "External safety" input operation	■ C 30
	■ Controlling the pump using communication software	■ C 41
	Maintenance	■ Precautions before maintenance
	■ Diagnosis and troubleshooting	■ D 20
	■ Pump running-in for ATH 31	■ D 30
	■ "Time Bearings" counter resetting.	■ D 40